



# **SureSelect Strand-Specific RNA Library Prep for Illumina Multiplexed Sequencing**

## **mRNA Library Preparation Protocol**

### **Protocol**

Version E0, March 2017

**SureSelect platform manufactured with Agilent  
SurePrint Technology**

**For Research Use Only. Not for use in diagnostic  
procedures.**



**Agilent Technologies**

## Notices

© Agilent Technologies, Inc. 2013-2015, 2017

No part of this manual may be reproduced in any form or by any means (including electronic storage and retrieval or translation into a foreign language) without prior agreement and written consent from Agilent Technologies, Inc. as governed by United States and international copyright laws.

### Manual Part Number

G9691-90010

### Edition

Version E0, March 2017

Printed in USA

Agilent Technologies, Inc.  
5301 Stevens Creek Blvd  
Santa Clara, CA 95051 USA

### Acknowledgement

Oligonucleotide sequences © 2006, 2008, and 2011 Illumina, Inc. All rights reserved. Only for use with the Illumina sequencer systems and associated assays.

### Technical Support

For technical product support, contact your local Agilent Support Services representative.

For US and Canada, call (800) 227-9770 (option 3,4,4). For other countries, find your support center telephone numbers at [www.agilent.com/chem/contactus](http://www.agilent.com/chem/contactus).

Or send an e-mail to:  
[SureSelect.Support@agilent.com](mailto:SureSelect.Support@agilent.com)

### Warranty

**The material contained in this document is provided “as is,” and is subject to being changed, without notice, in future editions. Further, to the maximum extent permitted by applicable law, Agilent disclaims all warranties, either express or implied, with regard to this manual and any information contained herein, including but not limited to the implied warranties of merchantability and fitness for a particular purpose. Agilent shall not be liable for errors or for incidental or consequential damages in connection with the furnishing, use, or performance of this document or of any information contained herein. Should Agilent and the user have a separate written agreement with warranty terms covering the material in this document that conflict with these terms, the warranty terms in the separate agreement shall control.**

### Technology Licenses

The hardware and/or software described in this document are furnished under a license and may be used or copied only in accordance with the terms of such license.

### Restricted Rights Legend

U.S. Government Restricted Rights. Software and technical data rights granted to the federal government include only those rights customarily provided to end user customers. Agilent provides this customary commercial license in Software and technical data pursuant to FAR 12.211 (Technical Data) and 12.212 (Computer Software) and, for the Department of Defense, DFARS 252.227-7015 (Technical Data - Commercial Items) and DFARS 227.7202-3 (Rights in Commercial Computer Software or Computer Software Documentation).

### Notice to Purchaser: Limited License

Use of this product is covered by one or more of the following US patents and corresponding patent claims outside the US: 6,258,569, 6,171,785, 6,127,155, 6,030,787, 5,994,056, 5,876,930, 5,804,375, 5,789,224, 5,773,258 (claims 1 and 6 only), 5,723,591, 5,677,152 (claims 1 to 23 only), 5,618,711, 5,538,848, and claims outside the US corresponding to expired US Patent No. 5,079,352. The purchase of this product includes a limited, non-transferable immunity from suit under the foregoing patent claims for using only this amount of product for the purchaser's own internal research. No right under any other patent claim and no right to perform commercial services of any kind, including without limitation reporting the results of purchaser's activities for a fee or other commercial consideration, is conveyed expressly, by implication, or by estoppel. This product is for research use only. Diagnostic uses under Roche patents require a separate license from Roche. Further information on purchasing licenses may be obtained by contacting the Director of Licensing, Applied Biosystems, 850 Lincoln Centre Drive, Foster City, California 94404, USA.

## Safety Notices

### CAUTION

A **CAUTION** notice denotes a hazard. It calls attention to an operating procedure, practice, or the like that, if not correctly performed or adhered to, could result in damage to the product or loss of important data. Do not proceed beyond a **CAUTION** notice until the indicated conditions are fully understood and met.

---

### WARNING

A **WARNING** notice denotes a hazard. It calls attention to an operating procedure, practice, or the like that, if not correctly performed or adhered to, could result in personal injury or death. Do not proceed beyond a **WARNING** notice until the indicated conditions are fully understood and met.

---

## In this Guide...

This guide describes an optimized protocol for Illumina paired-end multiplexed library preparation using the SureSelect Strand Specific RNA Library Prep system.

This protocol is specifically developed and optimized to prepare mRNA sequencing libraries from total RNA samples.

### **1 Before You Begin**

This chapter contains information (such as procedural notes, safety information, required reagents and equipment) that you should read and understand before you start an experiment.

### **2 Sample Preparation**

This chapter describes the steps to prepare cDNA sequencing libraries from mRNA isolated from total RNA samples.

### **3 Reference**

This chapter contains reference information, including component kit contents and index sequences.



## What's New in Version E0

- Updated storage duration and instructions for Actinomycin D stock solution (see [page 11](#), [page 17](#) and [page 22](#)).
- Support for sample processing in strip tubes (see [Table 2](#) on [page 12](#), with additional changes throughout protocol).
- Update to magnetic separator and low-adhesion microcentrifuge tubes supplier information (see [Table 2](#) on [page 12](#)).
- Updates to instructions for thawing kit components before use (see [page 17](#) and [page 28](#)).
- Updates to mixing instructions after addition of RNA Seq Bead Washing Buffer to sample mixtures (see [step 9](#) on [page 19](#) and [step 20](#) on [page 20](#)).
- Clarification to mixing instructions for kit-supplied master mixes, including vortexing duration and speed and pipetting considerations ([page 10](#), [page 22](#), [page 26](#), [page 28](#), [page 30](#), [page 32](#)).
- Updates to duration of certain bead collection steps (see [step 4](#) on [page 23](#) and [step 15](#) on [page 31](#)).
- Updates to instructions for bead drying and library elution during purification steps (for example, see [step 12](#) to [step 14](#) on [page 24](#)).
- Update to RNA fragmentation thermal cycler program to include 1-minute incubation at 4°C prior to indefinite hold at 4°C (see [Table 8](#) on [page 21](#)).
- Kit use-by information removed. See product labels and the Certificate of Analysis for each component for expiration date information.
- Updates to Bioanalyzer setup wording and link to DNA 1000 User Guide ([page 36](#)).
- Updates to supplier name for materials purchased from Thermo Fisher Scientific (see [Table 1](#) on [page 11](#)).
- Addition of [Optional Reagents and Equipment](#) table on [page 13](#).

- Updates to reagent descriptions to include vial format and cap color (for example, see [Table 4](#) on page 17).
- Addition of high-quality RNA control recommendation (see [Table 3](#) on page 13 and see [page 18](#)).
- Removal of reference information for obsolete kits containing indexing primers 1–96, provided in a clear plate (typically received before December 2014). If you need assistance with kits containing these obsolete indexing primer components, please contact [Technical Support](#).

# Content

<b>1</b>	<b>Before You Begin</b>	<b>9</b>
	Procedural Notes	10
	Safety Notes	10
	Required Reagents	11
	Required Equipment	12
	Optional Reagents and Equipment	13
<b>2</b>	<b>Sample Preparation</b>	<b>15</b>
	Step 1. Purify poly(A) RNA from total RNA	18
	Step 2. Fragment poly(A) RNA	21
	Step 3. Synthesize first-strand cDNA	22
	Step 4. Purify first strand cDNA using AMPure XP beads	24
	Step 5. Synthesize second-strand cDNA and repair the ends	26
	Step 6. Purify cDNA using AMPure XP beads	27
	Step 7. dA-Tail the cDNA 3' ends	28
	Step 8. Ligate adaptors	30
	Step 9. Purify adaptor-ligated DNA using AMPure XP beads	31
	Step 10. Amplify and index the adaptor-ligated cDNA library	32
	Step 11. Purify the amplified library with AMPure XP beads	35
	Step 12. Assess quality with the 2100 Bioanalyzer DNA 1000 Assay	36
	Step 13. Remove adaptor-dimers with AMPure XP beads (optional)	38
	Step 14. Prepare samples for multiplexed sequencing	40
	Step 15. Prepare and analyze sequencing samples	42
<b>3</b>	<b>Reference</b>	<b>45</b>
	Kit Contents	46
	Nucleotide Sequences of SureSelect <sup>XT</sup> Indexes A01 to H12	49

## Contents



# 1 Before You Begin

Procedural Notes	10
Safety Notes	10
Required Reagents	11
Optional Reagents and Equipment	13

Make sure you read and understand the information in this chapter and have the necessary equipment and reagents listed before you start an experiment.



## Procedural Notes

- To prevent contamination of reagents by nucleases, always wear powder-free laboratory gloves and use dedicated solutions and pipettors with nuclease-free aerosol-resistant tips.
- Maintain a clean work area.
- Avoid repeated freeze-thaw cycles of stock and diluted RNA and cDNA solutions. Possible stopping points, where samples may be stored at  $-20^{\circ}\text{C}$ , are marked in the protocol. Do not subject the samples to multiple freeze/thaw cycles.
- When preparing master mix reagent stock solutions for use:
  - 1 Thaw the reagent vial as rapidly as possible without heating above room temperature.
  - 2 Mix thoroughly on a vortex mixer at high speed for 5 seconds, then briefly spin in a centrifuge to drive the contents off of walls and lid.
  - 3 Store vials used during an experiment on ice or in a cold block.
  - 4 Library Preparation Master Mixes should not be frozen and thawed more than five times. If you plan to use the reagents in more than five experiments, aliquot to multiple vials to minimize freeze/thaw cycles for each vial.
- In general, follow Biosafety Level 1 (BL1) safety rules.

## Safety Notes

### CAUTION

- Wear appropriate personal protective equipment (PPE) when working in the laboratory.
-

## Required Reagents

**Table 1** Required Reagents for SureSelect Strand Specific RNA Library Prep

Description	Vendor and part number
SureSelect Strand Specific RNA Reagent Kit	Agilent
Illumina platforms (ILM), 16 Samples	p/n G9691A
Illumina platforms (ILM), 96 Samples	p/n G9691B
Actinomycin D*	Sigma p/n A1410
DMSO	Sigma p/n D8418
Agencourt AMPure XP Kit	Beckman Coulter Genomics
5 mL	p/n A63880
60 mL	p/n A63881
450 mL	p/n A63882
DNA 1000 Kit	Agilent p/n 5067-1504
1X Low TE Buffer (10 mM Tris-HCl, pH 8.0, 0.1 mM EDTA)	Thermo Fisher Scientific p/n 12090-015, or equivalent
100% Ethanol, molecular biology grade	Sigma-Aldrich p/n E7023

\* Actinomycin D should be obtained as a solid and prepared at 4  $\mu\text{g}/\mu\text{l}$  concentration in DMSO then stored in single-use aliquots at  $-20^{\circ}\text{C}$ , protected from light. The aliquots may be stored for up to one year before use. See [page 16](#) for additional information.

## Required Equipment

**Table 2** Required Equipment for SureSelect Strand Specific RNA Library Prep

Description	Vendor and part number
SureCycler 8800 Thermal Cycler, or equivalent	Agilent p/n G8800A
96 well plate module for SureCycler 8800 Thermal Cycler	Agilent p/n G8810A
SureCycler 8800-compatible plasticware:	
96-well plates	Agilent p/n 410088
OR	
8-well strip tubes	Agilent p/n 410092
Tube cap strips, domed	Agilent p/n 410096
2100 Bioanalyzer Laptop Bundle	Agilent p/n G2943CA
2100 Bioanalyzer Electrophoresis Set	Agilent p/n G2947CA
Centrifuge	Eppendorf Centrifuge model 5804 or equivalent
Low-Adhesion Tubes (RNase, DNase, and DNA-free)	USA Scientific
1.5 mL	p/n 1415-2600
0.5 mL	p/n 1405-2600
Magnetic separator	Thermo Fisher Scientific p/n 12331D or equivalent*
Multichannel pipette	Pipetman or equivalent
P10, P20, P200 and P1000 pipettes	Pipetman P10, P20, P200, P1000 or equivalent
Vortex mixer	
Ice bucket	
Powder-free gloves	
Sterile, nuclease-free aerosol barrier pipette tips	

\* Select a magnetic separator compatible with the 96-well plates or 8-well strip tubes used for sample processing and configured to collect magnetic particles on one side of each well. Do not use a magnetic separator configured to collect the particles in a ring formation.




## Optional Reagents and Equipment

**Table 3** Optional Reagents and Equipment

Description	Vendor and part number
Labnet MPS1000 Mini Plate Spinner	Labnet International p/n C1000
Universal Human Reference RNA	Agilent p/n 740000

**1 Before You Begin**  
Optional Reagents and Equipment

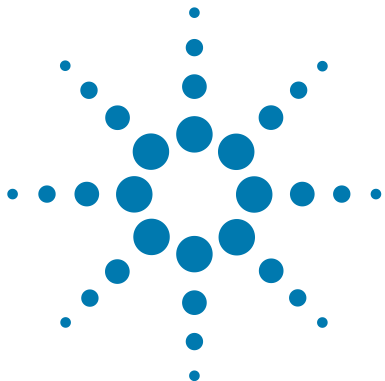

## 2 Sample Preparation


- Step 1. Purify poly(A) RNA from total RNA 18
- Step 2. Fragment poly(A) RNA 21
- Step 3. Synthesize first-strand cDNA 22
- Step 4. Purify first strand cDNA using AMPure XP beads 24
- Step 5. Synthesize second-strand cDNA and repair the ends 26
- Step 6. Purify cDNA using AMPure XP beads 27
- Step 7. dA-Tail the cDNA 3' ends 28
- Step 8. Ligate adaptors 30
- Step 9. Purify adaptor-ligated DNA using AMPure XP beads 31
- Step 10. Amplify and index the adaptor-ligated cDNA library 32
- Step 11. Purify the amplified library with AMPure XP beads 35
- Step 12. Assess quality with the 2100 Bioanalyzer DNA 1000 Assay 36
- Step 13. Remove adaptor-dimers with AMPure XP beads (optional) 38
- Step 14. Prepare samples for multiplexed sequencing 40
- Step 15. Prepare and analyze sequencing samples 42

See [Figure 1](#) for a summary of the SureSelect mRNA sequencing workflow.

This section contains instructions for strand-specific RNA sequencing cDNA library preparation for the Illumina platform.



# SureSelect Strand-Specific mRNA Library Preparation for NGS Workflow

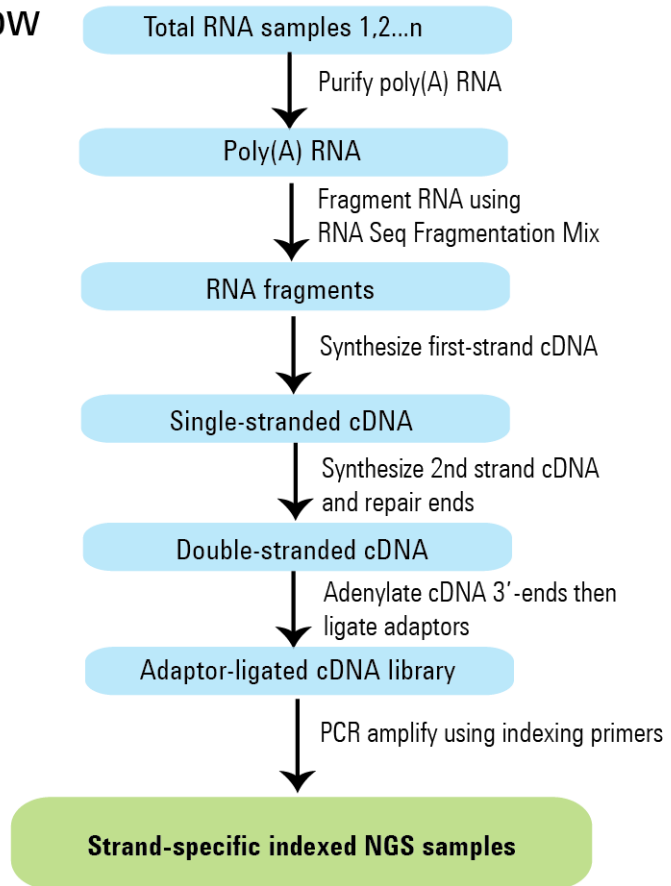

**Figure 1** Overall mRNA sequencing sample preparation workflow.

Before you begin, prepare a stock solution of 4 µg/µL Actinomycin D in DMSO. Aliquot the stock solution into single-use volumes (typically 3 µL). Store the aliquots at –20°C, protected from light. Do not subject the aliquots to multiple freeze-thaw cycles. The aliquots may be stored for up to one year before use in the library preparation protocol. During the library preparation protocol, an aliquot of the DMSO stock is diluted with water, immediately before use, to a final Actinomycin D concentration of 120 ng/µL. (See [page 22](#) for more information.)

Just before starting the protocol, bring the reagents listed in [Table 4](#) to room temperature and thaw the reagents listed in [Table 5](#) on ice.

**Table 4** Reagents brought to room temperature before use in protocol

Kit Component	Storage Location	Where Used in Protocol
Oligo(dT) Microparticles (tube with brown cap or bottle)	RNA Library Prep Kit Box 2, 4°C	<a href="#">page 18</a>
RNA Seq Bead Washing Buffer (bottle)	RNA Library Prep Kit Box 2, 4°C	<a href="#">page 19</a>
RNA Seq Bead Elution Buffer (tube with green cap or bottle)	RNA Library Prep Kit Box 2, 4°C	<a href="#">page 20</a>
RNA Seq Bead Binding Buffer (tube with purple cap or bottle)	RNA Library Prep Kit Box 2, 4°C	<a href="#">page 20</a>
Nuclease Free Water (bottle)	RNA Library Prep Kit Box 2, 4°C	<a href="#">page 18</a>

**Table 5** Reagents thawed and held on ice before use in protocol

Kit Component	Storage Location	Where Used in Protocol
RNA Seq Fragmentation Mix (tube with red cap or bottle)	RNA Library Prep Kit Box 1, –20°C	<a href="#">page 21</a>
RNA Seq First Strand Master Mix (tube with orange cap)	RNA Library Prep Kit Box 1, –20°C	<a href="#">page 22</a>
RNA Seq Second-Strand + End-Repair Enzyme Mix (tube with blue cap or bottle)	RNA Library Prep Kit Box 1, –20°C	<a href="#">page 26</a>
RNA Seq Second-Strand + End-Repair Oligo Mix (tube with yellow cap)	RNA Library Prep Kit Box 1, –20°C	<a href="#">page 26</a>

## 2 Sample Preparation

### Step 1. Purify poly(A) RNA from total RNA

#### Step 1. Purify poly(A) RNA from total RNA

In this step, poly(A) RNA is purified from total RNA using two serial rounds of binding to oligo(dT) magnetic particles.

Before you begin, prepare total RNA for each sample in the run. The amount of total RNA needed for the library preparation protocol is 50 ng to 4 µg.

#### NOTE

For optimal performance, total RNA samples should have an RNA Integrity Number (RIN) of 8 or more, based on analysis using Agilent's 2100 Bioanalyzer.

Consider preparing an additional sequencing library in parallel, using a high-quality control RNA sample, such as Agilent's Universal Human Reference RNA (p/n 740000). Use of this control is especially recommended during the first run of the protocol, to verify that all protocol steps are being successfully performed. Routine use of this control is helpful for any required troubleshooting, in order to differentiate any performance issues related to RNA input from other factors.

Make sure the reagents to be used in the protocol have been thawed and kept at the appropriate temperature, as indicated in [Table 4](#) and [Table 5](#) on page 17.

- 1 Prepare each total RNA sample in a final volume of 25 µL of nuclease-free water and place the samples in separate wells of a 96-well plate or strip tube.
- 2 Vortex the Oligo(dT) Microparticles until the suspension appears homogeneous and consistent in color. If bead aggregates are still present after vortexing, mix thoroughly by pipetting up and down until the suspension appears homogeneous.
- 3 Add 25 µL of the homogeneous Oligo(dT) bead suspension to each total RNA sample well.
- 4 Seal the wells, then gently vortex for 5 seconds and briefly spin the plate or strip tube in a centrifuge or mini-plate spinner to collect the liquid without pelleting the beads.

- 5** Incubate the plate or strip tube in the SureCycler thermal cycler (with the heated lid ON) and run the program in [Table 6](#) to denature the RNA.

**Table 6** Thermal cycler program for RNA denaturation

Step	Temperature	Time
Step 1	65°C	5 minutes
Step 2	4°C	1 minute
Step 3	4°C	Hold

- 6** After the thermal cycler reaches the 4°C Hold step, remove the plate or strip tube and incubate at room temperature for 5 minutes, to allow poly(A) RNA binding to the oligo(dT) beads.
- 7** Move the plate or strip tube to a magnetic separation device at room temperature. Wait for the solution to clear (approximately 2 to 5 minutes).
- 8** While keeping the plate or strip tube in the magnetic stand, carefully remove and discard the cleared solution from each well. Do not touch or disturb the beads while removing the solution.
- 9** Remove the plate or strip tube from the magnetic stand. Gently add 200 µL of RNA Seq Bead Washing Buffer to each well.  
Mix by pipetting up and down 10 times, using a P200 pipette set to 150 µL, without introducing bubbles.

### CAUTION

The Bead Washing Buffer contains detergent. It is important to process mixtures of the beads with the wash buffer without introducing bubbles or foam. If bubbles or foam are present during the wash steps, briefly spin the plate or strip tube in a centrifuge before continuing.

- 10** Place the plate or strip tube in the magnetic stand at room temperature. Wait for the solution to clear (at least 2 minutes).
- 11** While keeping the plate or strip tube in the magnetic stand, carefully remove and discard the cleared solution from each well. Do not touch or disturb the beads while removing the solution.

## 2 Sample Preparation

### Step 1. Purify poly(A) RNA from total RNA

- 12** Remove the plate or strip tube from the magnetic stand. Add 25  $\mu$ L of RNA Seq Bead Elution Buffer to each sample well.
- 13** Seal the wells, then gently vortex for 5 seconds. Briefly spin the plate or strip tube in a centrifuge or mini-plate spinner to collect the liquid.
- 14** Incubate the plate or strip tube in the thermal cycler (with the heated lid ON) and run the program in [Table 7](#).

**Table 7** Thermal cycler program for RNA elution

Step	Temperature	Time
Step 1	80°C	2 minutes
Step 2	4°C	1 minute
Step 3	4°C	Hold

- 15** After the thermal cycler reaches the 4°C Hold step, remove the plate or strip tube and add 25  $\mu$ L of RNA Seq Bead Binding Buffer to each sample well.
- 16** Seal the wells, then gently vortex for 5 seconds. Briefly spin the plate or strip tube in a centrifuge or mini-plate spinner to collect the liquid.
- 17** Incubate the samples at room temperature for 5 minutes, to allow poly(A) RNA to re-bind the beads.
- 18** Place the plate or strip tube in the magnetic stand at room temperature. Wait for the solution to clear (at least 2 minutes).
- 19** While keeping the plate or strip tube in the magnetic stand, carefully remove and discard the cleared solution from each well. Do not touch or disturb the beads while removing the solution.
- 20** Remove the plate or strip tube from the magnetic stand. Gently add 200  $\mu$ L of RNA Seq Bead Washing Buffer to each well.  
Mix by pipetting up and down 10 times, using a P200 pipette set to 150  $\mu$ L, without introducing bubbles. If bubbles or foam are present, spin the plate or strip tube briefly before continuing.
- 21** Place the plate or strip tube in the magnetic stand at room temperature. Wait for the solution to clear (at least 2 minutes).
- 22** While keeping the plate or strip tube in the magnetic stand, carefully remove and discard the cleared solution from each well. Do not touch or disturb the beads while removing the solution.
- 23** Proceed immediately to [Step 2. Fragment poly\(A\) RNA](#).



## Step 2. Fragment poly(A) RNA

In this step, the purified poly(A) RNA is chemically-fragmented to a size appropriate for RNA sequencing library preparation.

- 1** Remove the plate or strip tube, containing the collected poly(A) RNA-bound beads, from the magnetic stand.
- 2** To each sample well add 19  $\mu$ L of RNA Seq Fragmentation Mix.
- 3** Seal the wells, then gently vortex for 5 seconds. Briefly spin the plate or strip tube in a centrifuge or mini-plate spinner to collect the liquid.
- 4** Incubate the plate or strip tube in the thermal cycler (with the heated lid ON) and run the program in [Table 8](#).

**Table 8** Thermal cycler program for RNA fragmentation

Step	Temperature	Time
Step 1	94°C	8 minutes
Step 2	4°C	1 minute
Step 3	4°C	Hold

- 5** During the 8-minute incubation at 94°C, complete the first-strand cDNA reagent preparation steps ([step 1](#) through [step 3](#)) on [page 22](#).

## Step 3. Synthesize first-strand cDNA

### CAUTION

To ensure strand-specificity, you must prepare the 120 ng/ $\mu$ L Actinomycin D solution in [step 1](#), below, immediately before use. The stock solution of 4  $\mu$ g/ $\mu$ L Actinomycin D in DMSO must be prepared less than one year prior to use and stored in single-use aliquots at  $-20^{\circ}\text{C}$ , protected from light.

- 1 Prepare a fresh 120 ng/ $\mu$ L Actinomycin D dilution in water, using an aliquot of the 4  $\mu$ g/ $\mu$ L Actinomycin D stock solution in DMSO, according to [Table 9](#).

**Table 9** Preparation of 120 ng/ $\mu$ L Actinomycin D

Reagent	Volume for up to 96-reaction run (includes excess)
Actinomycin D (4 $\mu$ g/ $\mu$ L in DMSO)	3 $\mu$ L
Nuclease-free water	97 $\mu$ L
<b>Total</b>	<b>100 <math>\mu</math>L</b>

### CAUTION

The RNA Seq First Strand Master Mix used in this step is viscous. Mix thoroughly by vortexing at high speed for 5 seconds before removing an aliquot for use and after combining with other solutions. Pipetting up and down is not sufficient to mix this reagent.

- 2 Vortex the thawed vial of RNA Seq First Strand Master Mix 5 seconds at high speed to ensure homogeneity.

- 3 Prepare the appropriate amount of RNA Seq First Strand Master Mix + Actinomycin D mixture according to [Table 10](#) below. Mix by vortexing, then spin briefly and keep on ice.

**Table 10** Preparation of First Strand Master Mix/Actinomycin D mixture

Reagent	Volume for 1 reaction	Volume for 16 reactions (includes excess)
Actinomycin D (120 ng/μl in H <sub>2</sub> O)	0.5 μL	8.5 μL
RNA Seq First Strand Master Mix (orange cap)	8.0 μL	136 μL
<b>Total</b>	<b>8.5 μL</b>	<b>144.5 μL</b>

- 4 Once the RNA fragmentation thermal cycler program has reached the 4°C Hold step (Step 3 in [Table 8](#)), transfer the sample plate or strip tube from the thermal cycler to the magnetic stand at room temperature. Leave the plate or strip tube on the magnetic stand for at least 2 minutes, until the solution is clear.
- 5 Keep the RNA sample plate or strip tube in the magnetic stand at room temperature while you carefully transfer 17 μL of each supernatant to a fresh well. Do not touch or disturb the beads while removing the fragmented RNA solution. Once the supernatant from all wells has been transferred, place the samples on ice or in a cold block.
- 6 Add 8.5 μL of First Strand Master Mix/Actinomycin D mixture prepared in [step 2](#) to each RNA sample well.
- 7 Seal the wells. Mix thoroughly by vortexing the plate or strip tube for 5 seconds or by agitating the plate or strip tube on a plate shaker at approximately 1600 rpm for 5 seconds.
- 8 Spin the plate or strip tube briefly to collect the liquid.
- 9 Incubate the plate or strip tube in the thermal cycler (with the heated lid ON) and run the program in [Table 11](#).

**Table 11** Thermal cycler program for first-strand cDNA synthesis

Step	Temperature	Time
Step 1	25°C	10 minutes
Step 2	37°C	40 minutes
Step 3	4°C	Hold

## 2 Sample Preparation

### Step 4. Purify first strand cDNA using AMPure XP beads

#### Step 4. Purify first strand cDNA using AMPure XP beads

- 1 Let the AMPure XP beads come to room temperature for at least 30 minutes. *Do not freeze the beads at any time.*
- 2 Prepare 400  $\mu\text{L}$  of ~~SURE~~ 70% ethanol per sample, plus excess, for use in step 8.

#### NOTE

The freshly-prepared 70% ethanol may be used for subsequent purification steps run on the same day. The complete SureSelect RNA Library Preparation protocol requires 1.6 mL of fresh 70% ethanol per sample.

- 3 Mix the bead suspension well so that the suspension appears homogeneous and consistent in color.
- 4 Add 46  $\mu\text{L}$  of the homogeneous bead suspension to each 25.5- $\mu\text{L}$  sample in the PCR plate or strip tube. Seal the wells, then vortex the plate or strip tube for 5 seconds. Briefly spin to collect the liquid.
- 5 Incubate samples for 5 minutes at room temperature.
- 6 Put the plate or strip tube on the magnetic stand at room temperature. Wait for the solution to clear (at least 5 minutes).
- 7 While keeping the plate or strip tube in the magnetic stand, carefully remove and discard the cleared solution from each well. Do not touch the beads while removing the solution.
- 8 Continue to keep the plate or strip tube in the magnetic stand while you dispense 200  $\mu\text{L}$  of fresh 70% ethanol in each sample well.
- 9 Wait for 1 minute to allow any disturbed beads to settle, then remove the ethanol.
- 10 Repeat [step 8](#) and [step 9](#) step once for a total of two washes.
- 11 After removing the 200  $\mu\text{L}$  ethanol supernatant from the second wash, spin the plate or strip tube briefly, return the samples to the magnetic stand, and then remove any remaining ethanol droplets with a pipette.
- 12 Dry the beads on the thermal cycler (with lid open) at 37°C for up to 1 minute. Do not overdry the beads.
- 13 Add 21  $\mu\text{L}$  nuclease-free water to each sample well.
- 14 Seal the wells, then vortex the plate or strip tube for 5 seconds. Verify that all beads have been resuspended, with no visible clumps in the suspension or bead pellets retained on the sides of the wells. Briefly spin to collect the liquid.
- 15 Incubate for 2 minutes at room temperature.

## Step 4. Purify first strand cDNA using AMPure XP beads

- 16** Put the plate or strip tube in the magnetic stand and leave for 2 minutes or until the solution is clear.
- 17** Remove 20  $\mu$ L of cleared supernatant to a fresh well and keep on ice. You can discard the beads at this time.
- 18** Proceed immediately to “[Step 5. Synthesize second-strand cDNA and repair the ends](#)” on page 26.

## 2 Sample Preparation

### Step 5. Synthesize second-strand cDNA and repair the ends

## Step 5. Synthesize second-strand cDNA and repair the ends

#### CAUTION

The RNA Seq Second Strand + End Repair Enzyme Mix used in this step is viscous. Mix thoroughly by vortexing at high speed for 5 seconds before removing an aliquot for use and after combining with other solutions. Pipetting up and down is not sufficient to mix this reagent.

- 1 Vortex the thawed vials of RNA Seq Second Strand + End Repair Enzyme Mix and RNA Seq Second Strand + End Repair Oligo Mix at high speed for 5 seconds to ensure homogeneity.
- 2 Add 25  $\mu\text{L}$  of RNA Seq Second Strand + End Repair Enzyme Mix to each 20- $\mu\text{L}$  purified first-strand cDNA sample. Keep on ice.
- 3 Add 5  $\mu\text{L}$  of RNA Seq Second Strand + End Repair Oligo Mix to each sample well, for a total reaction volume of 50  $\mu\text{L}$ . Keep on ice.
- 4 Seal the wells. Mix thoroughly by vortexing the plate or strip tube for 5 seconds or by agitating the plate or strip tube on a plate shaker at approximately 1600 rpm for 5 seconds.
- 5 Spin the plate or strip tube briefly to collect the liquid.
- 6 Incubate the plate or strip tube in the thermal cycler and run the program in [Table 12](#). **Do not use a heated lid.**

**Table 12** Thermal cycler program for second-strand synthesis and end repair

Step	Temperature	Time
Step 1	16°C	30 minutes
Step 2	4°C	Hold

## Step 6. Purify cDNA using AMPure XP beads

- 1 Let the AMPure XP beads come to room temperature for at least 30 minutes. *Do not freeze the beads at any time.*
- 2 Prepare 400  $\mu\text{L}$  of fresh 70% ethanol per sample, plus excess, for use in [step 8](#).
- 3 Mix the bead suspension well so that the suspension appears homogeneous and consistent in color.
- 4 Add 90  $\mu\text{L}$  of the homogeneous bead suspension to each 50- $\mu\text{L}$  sample in the PCR plate or strip tube. Seal the wells, then vortex the plate or strip tube for 5 seconds. Briefly spin to collect the liquid.
- 5 Incubate samples for 5 minutes at room temperature.
- 6 Put the plate or strip tube on the magnetic stand at room temperature. Wait for the solution to clear (at least 5 minutes).
- 7 While keeping the plate or strip tube in the magnetic stand, carefully remove and discard the cleared solution from each well. Do not touch the beads while removing the solution.
- 8 Continue to keep the plate or strip tube in the magnetic stand while you dispense 200  $\mu\text{L}$  of fresh 70% ethanol in each sample well.
- 9 Wait for 1 minute to allow any disturbed beads to settle, then remove the ethanol.
- 10 Repeat [step 8](#) and [step 9](#) step once for a total of two washes.
- 11 Dry the beads on the thermal cycler (with lid open) at 37°C for up to 2 minutes. Do not overdry the beads.
- 12 Add 21  $\mu\text{L}$  nuclease-free water to each sample well.
- 13 Seal the wells, then vortex the plate or strip tube for 5 seconds. Verify that all beads have been resuspended, with no visible clumps in the suspension or bead pellets retained on the sides of the wells. Briefly spin to collect the liquid.
- 14 Incubate for 2 minutes at room temperature.
- 15 Put the plate or strip tube in the magnetic stand and leave for 2 minutes or until the solution is clear.
- 16 Remove 20  $\mu\text{L}$  of cleared supernatant to a fresh well and keep the samples on ice. You can discard the beads at this time.

**Stopping Point** If you do not continue to the next step, seal the wells and store at  $-20^{\circ}\text{C}$ .

## Step 7. dA-Tail the cDNA 3' ends

Before continuing the protocol, thaw reagents listed in [Table 13](#) (used for the remainder of library preparation steps) and keep on ice.

**Table 13** Reagents thawed and held on ice before use in protocol

Kit Component	Storage Location	Where Used in Protocol
RNA Seq dA Tailing Master Mix (tube with green cap or bottle)	RNA Library Prep Kit Box 1, –20°C	<a href="#">page 28</a>
SureSelect Ligation Master Mix (tube with purple cap)	RNA Library Prep Kit Box 1, –20°C	<a href="#">page 30</a>
SureSelect Oligo Adaptor Mix (tube with blue cap)	RNA Library Prep Kit Box 1, –20°C	<a href="#">page 30</a>
RNA Seq ILM Reverse PCR Primer (tube with black cap)	RNA Library Prep Kit Box 1, –20°C	<a href="#">page 32</a>
RNA Seq PCR Master Mix (tube with red cap or bottle)	RNA Library Prep Kit Box 1, –20°C	<a href="#">page 32</a>
Uracil DNA Glycosylase (UDG) (tube with yellow cap)	RNA Library Prep Kit Box 1, –20°C	<a href="#">page 33</a>
SureSelect Primer (tube with brown cap)	RNA Library Prep Kit Box 1, –20°C	<a href="#">page 33</a>
SureSelect <sup>XT</sup> Indexes, 8 bp (tubes with white caps or blue 96-well plate)	RNA Library Prep Kit Box 1, –20°C	<a href="#">page 33</a>

### CAUTION

The RNA Seq dA Tailing Master Mix used in this step is viscous. Mix thoroughly by vortexing at high speed for 5 seconds before removing an aliquot for use and after combining with other solutions. Pipetting up and down is not sufficient to mix this reagent.

- 1 Vortex the thawed vial of RNA Seq dA Tailing Master Mix 5 seconds at high speed to ensure homogeneity.
- 2 Add 20 µL of RNA Seq dA Tailing Master Mix to each 20-µL purified, end-repaired cDNA sample.



Pipette the master mix slowly to ensure that the full volume is dispensed.

- 3 Seal the wells. Mix thoroughly by vortexing the plate or strip tube for 5 seconds or by agitating the plate or strip tube on a plate shaker at approximately 1600 rpm for 5 seconds.
- 4 Spin the plate or strip tube briefly to collect the liquid.
- 5 Incubate the plate or strip tube in the thermal cycler and run the program in [Table 14](#). **Do not use a heated lid.**

**Table 14** Thermal cycler program for dA-tailing

Step	Temperature	Time
Step 1	37°C	15 minutes
Step 2	4°C	Hold

## Step 8. Ligate adaptors

### CAUTION

The SureSelect Ligation Master Mix used in this step is viscous. Mix thoroughly by vortexing at high speed for 5 seconds before removing an aliquot for use and after combining with other solutions. Pipetting up and down is not sufficient to mix this reagent.

- 1 Vortex the thawed vials of SureSelect Ligation Master Mix and SureSelect Oligo Adaptor Mix at high speed for 5 seconds to ensure homogeneity.
- 2 Transfer the cDNA samples to ice, then add 5  $\mu$ L of SureSelect Ligation Master Mix to each A-tailed cDNA sample well.  
Pipette the master mix slowly to ensure that the full volume is dispensed.
- 3 Add 5  $\mu$ L of SureSelect Oligo Adaptor Mix to each sample.
- 4 Seal the wells. Mix thoroughly by vortexing the plate or strip tube for 5 seconds or by agitating the plate or strip tube on a plate shaker at approximately 1600 rpm for 5 seconds.
- 5 Spin the plate or strip tube briefly to collect the liquid.
- 6 Incubate the plate or strip tube in the thermal cycler and run the program in [Table 15](#). **Do not use a heated lid.**

**Table 15** Thermal cycler program for adaptor ligation

Step	Temperature	Time
Step 1	20°C	15 minutes
Step 2	4°C	Hold

## Step 9. Purify adaptor-ligated DNA using AMPure XP beads

- 1 Let the AMPure XP beads come to room temperature for at least 30 minutes. *Do not freeze the beads at any time.*
- 2 Prepare 400  $\mu\text{L}$  of fresh 70% ethanol per sample, plus excess, for use in [step 8](#).
- 3 Mix the bead suspension well so that the suspension appears homogeneous and consistent in color.
- 4 Add 60  $\mu\text{L}$  of the homogeneous bead suspension to each 50- $\mu\text{L}$  sample in the PCR plate or strip tube. Seal the wells, then vortex the plate or strip tube for 5 seconds. Briefly spin to collect the liquid.
- 5 Incubate samples for 5 minutes at room temperature.
- 6 Put the plate or strip tube on the magnetic stand at room temperature. Wait for the solution to clear (at least 5 minutes).
- 7 While keeping the plate or strip tube in the magnetic stand, carefully remove and discard the cleared solution from each well. Do not touch the beads while removing the solution.
- 8 Continue to keep the plate or strip tube in the magnetic stand while you dispense 200  $\mu\text{L}$  of fresh 70% ethanol in each sample well.
- 9 Wait for 1 minute to allow any disturbed beads to settle, then remove the ethanol.
- 10 Repeat [step 8](#) and [step 9](#) step once for a total of two washes.
- 11 Dry the beads on the thermal cycler (with lid open) at 37°C for up to 3 minutes. Do not overdry the beads.
- 12 Add 18  $\mu\text{L}$  nuclease-free water to each sample well.
- 13 Seal the wells, then vortex the plate or strip tube for 5 seconds. Verify that all beads have been resuspended, with no visible clumps in the suspension or bead pellets retained on the sides of the wells. Briefly spin to collect the liquid.
- 14 Incubate for 2 minutes at room temperature.
- 15 Put the plate or strip tube in the magnetic stand and leave for 2 minutes or until the solution is clear.
- 16 Remove 17  $\mu\text{L}$  of cleared supernatant to a fresh well and keep on ice. You can discard the beads at this time.

## 2 Sample Preparation

### Step 10. Amplify and index the adaptor-ligated cDNA library

## Step 10. Amplify and index the adaptor-ligated cDNA library

In this step, the adaptor ligated cDNA is amplified in a three-primer PCR that includes the appropriate indexing primer. Amplification cycle number is based on the initial amount of RNA sample used for library preparation (see [Table 19](#) on page 34).

#### NOTE

The PCR protocol detailed below is appropriate for mRNA library amplification and PCR indexing without target enrichment. If samples will be target-enriched after library preparation, see publication part number G9691-90000 at [www.genomics.agilent.com](http://www.genomics.agilent.com) for the appropriate pre-capture amplification and post-capture indexing protocols.

- 1 Prepare a 1:20 dilution of the RNA Seq ILM Reverse PCR Primer, according to [Table 16](#).

**Table 16** Preparation of reverse primer dilution

Reagent	Volume for up to 16 reactions (includes excess)
Nuclease-free water	19 $\mu$ L
RNA Seq ILM Reverse PCR Primer (tube with black cap)	1 $\mu$ L
<b>Total</b>	<b>20 <math>\mu</math>L</b>

#### CAUTION

The RNA Seq PCR Master Mix used at this step is highly viscous and thorough mixing is critical for optimal kit performance. Mix by vortexing at high speed for 5 seconds before removing an aliquot for use and after combining with other solutions in [Table 17](#). Pipetting up and down is not sufficient to mix this reagent.

- 2 Vortex the thawed vial of RNA Seq PCR Master Mix 5 seconds at high speed to ensure homogeneity.

## Step 10. Amplify and index the adaptor-ligated cDNA library

- 3** Prepare the appropriate volume of PCR reaction mix, as described in [Table 17](#), on ice. Mix well by vortexing at high speed then spin briefly to collect the liquid.

**Table 17** Preparation of PCR Reaction Mix

Reagent	Volume for 1 reaction	Volume for 16 reactions (includes excess)
RNA Seq PCR Master Mix (tube with red cap or bottle)	25 $\mu$ L	425 $\mu$ L
Uracil DNA Glycosylase (UDG) (tube with yellow cap)	1 $\mu$ L	17 $\mu$ L
SureSelect Primer (forward primer; tube with brown cap)	1 $\mu$ L	17 $\mu$ L
1:20 dilution of RNA Seq ILM Reverse PCR Primer prepared in <a href="#">Table 16</a>	1 $\mu$ L	17 $\mu$ L
<b>Total</b>	<b>28 <math>\mu</math>L</b>	<b>476 <math>\mu</math>L</b>

- 4** Add 28  $\mu$ L of the PCR reaction mix prepared in [step](#) to each sample well containing 17  $\mu$ L of purified, adaptor-ligated cDNA. Mix by pipetting. Change pipette tips between samples.
- 5** Add 5  $\mu$ L of the appropriate indexing primer (Index A01–H06) to each PCR reaction mixture well.

**CAUTION**

For 96-reaction kits, only use the indexing primers provided in columns 1 through 6 (Indexes A01–H06) in [step 5](#). Provided plates also contain indexing primers in columns 7 through 12, which are used in other SureSelect protocols.

See [page 48](#) for a plate map.

- 6** Seal the wells. Mix thoroughly by vortexing the plate or strip tube for 5 seconds or by agitating the plate or strip tube on a plate shaker at approximately 1600 rpm for 5 seconds.
- 7** Spin the plate or strip tube briefly to collect the liquid.

## 2 Sample Preparation

### Step 10. Amplify and index the adaptor-ligated cDNA library

- 8 Incubate the plate or strip tube in the thermal cycler (with the heated lid ON) and run the program in [Table 18](#).

**Table 18** Thermal cycler program for mRNA Library PCR indexing

Segment	Number of Cycles	Temperature	Time
1	1	37°C	15 minutes
2	1	95°C	2 minutes
3	10–16 cycles (see <a href="#">Table 19</a> )	95°C	30 seconds
		65°C	30 seconds
		72°C	1 minute
4	1	72°C	5 minutes
5	1	4°C	Hold

**Table 19** mRNA Library PCR indexing cycle number recommendations

Amount of total RNA used for library prep	Cycle Number
50 ng–200 ng	14–16
201 ng–2 µg	12–14
2.1 µg–4 µg	10–12

#### NOTE

If you started with the minimum total RNA input amount of 50 ng, use 16 amplification cycles in the PCR indexing amplification program ([Table 18](#)).

## Step 11. Purify the amplified library with AMPure XP beads

- 1 Let the AMPure XP beads come to room temperature for at least 30 minutes. *Do not freeze the beads at any time.*
- 2 Prepare 400  $\mu\text{L}$  of fresh 70% ethanol per sample, plus excess, for use in [step 8](#).
- 3 Mix the bead suspension well so that the suspension appears homogeneous and consistent in color.
- 4 Add 60  $\mu\text{L}$  of the homogeneous bead suspension to each 50- $\mu\text{L}$  PCR reaction in the PCR plate or strip tube. Seal the wells, then vortex the plate or strip tube for 5 seconds. Briefly spin to collect the liquid.
- 5 Incubate samples for 5 minutes at room temperature.
- 6 Put the plate or strip tube on the magnetic stand at room temperature. Wait for the solution to clear (at least 5 minutes).
- 7 While keeping the plate or strip tube in the magnetic stand, carefully remove and discard the cleared solution from each well. Do not touch the beads while removing the solution.
- 8 Continue to keep the plate or strip tube in the magnetic stand while you dispense 200  $\mu\text{L}$  of fresh 70% ethanol in each sample well.
- 9 Wait for 1 minute to allow any disturbed beads to settle, then remove the ethanol.
- 10 Repeat [step 8](#) and [step 9](#) step once for a total of two washes.
- 11 Dry the beads on the thermal cycler (with lid open) at 37°C for up to 3 minutes. Do not overdry the beads.
- 12 Add 26  $\mu\text{L}$  nuclease-free water to each sample well.
- 13 Seal the wells, then vortex the plate or strip tube for 5 seconds. Verify that all beads have been resuspended, with no visible clumps in the suspension or bead pellets retained on the sides of the wells. Briefly spin to collect the liquid.
- 14 Incubate for 2 minutes at room temperature.
- 15 Put the plate or strip tube in the magnetic stand and leave for 2 minutes or until the solution is clear.
- 16 Remove 25  $\mu\text{L}$  of cleared supernatant to a fresh well and keep on ice. You can discard the beads at this time.

**Stopping Point** If you do not continue to the next step, seal the wells and store at  $-20^{\circ}\text{C}$ .

## Step 12. Assess quality with the 2100 Bioanalyzer DNA 1000 Assay

Use a Bioanalyzer DNA 1000 chip and reagent kit. See the *DNA 1000 Kit Guide* at [www.genomics.agilent.com](http://www.genomics.agilent.com) for more information on doing this step.


- 1 Check that the 2100 Bioanalyzer electrodes have been cleaned as instructed in the reagent kit guide.
- 2 Open the 2100 Expert Software (version B.02.02 or higher), turn on the 2100 Bioanalyzer and check communication.
- 3 Prepare the chip, samples and ladder as instructed in the reagent kit guide, using 1  $\mu$ L of each sample for the analysis.
- 4 Load the prepared chip into the 2100 Bioanalyzer. Within the instrument context, choose the DNA 1000 assay from the drop down list. Start the run within five minutes after preparation.
- 5 Verify the results. A sample electropherogram is shown in [Figure 2](#).
  - a Measure the concentration of the library by integrating under the peak at approximately 200 to 600 bp. For accurate quantification, make sure that the concentration falls within the linear range of the assay.
  - b Analyze the 100 to 150-bp range of the electropherogram. A peak at this position indicates the presence of adaptor-dimers in the cDNA sample, which need to be removed for optimal sequencing results.

If adaptor-dimers are observed, dilute the prepared library sample to a final volume of 50  $\mu$ L using nuclease-free water, then proceed to “[Step 13. Remove adaptor-dimers with AMPure XP beads \(optional\)](#)” on page 38.

If adaptor-dimers are not observed, proceed to “[Step 14. Prepare samples for multiplexed sequencing](#)” on page 40.



## Step 12. Assess quality with the 2100 Bioanalyzer DNA 1000 Assay

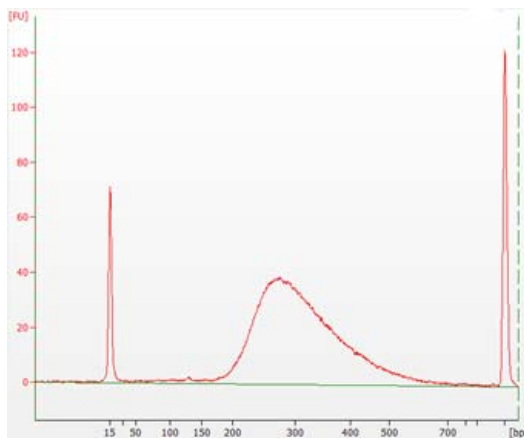

**Figure 2** Analysis of amplified prepped library cDNA using a DNA 1000 assay. The electropherogram shows a single peak in the size range of 200 to 600 bp.

## 2 Sample Preparation

### Step 13. Remove adaptor-dimers with AMPure XP beads (optional)

#### Step 13. Remove adaptor-dimers with AMPure XP beads (optional)

- 1 Let the AMPure XP beads come to room temperature for at least 30 minutes. *Do not freeze the beads at any time.*
- 2 Prepare 400  $\mu\text{L}$  of fresh 70% ethanol per sample, plus excess, for use in [step 8](#).
- 3 Mix the bead suspension well so that the suspension appears homogeneous and consistent in color.
- 4 Add 60  $\mu\text{L}$  of the homogeneous bead suspension to each 50- $\mu\text{L}$  library sample in the PCR plate or strip tube. Seal the wells, then vortex the plate or strip tube for 5 seconds. Briefly spin to collect the liquid.
- 5 Incubate samples for 5 minutes at room temperature.
- 6 Put the plate or strip tube on the magnetic stand at room temperature. Wait for the solution to clear (at least 5 minutes).
- 7 While keeping the plate or strip tube in the magnetic stand, carefully remove and discard the cleared solution from each well. Do not touch the beads while removing the solution.
- 8 Continue to keep the plate or strip tube in the magnetic stand while you dispense 200  $\mu\text{L}$  of fresh 70% ethanol in each sample well.
- 9 Wait for 1 minute to allow any disturbed beads to settle, then remove the ethanol.
- 10 Repeat [step 8](#) and [step 9](#) step once for a total of two washes.
- 11 Dry the beads on the thermal cycler (with lid open) at 37°C for up to 3 minutes. Do not overdry the beads.
- 12 Add 26  $\mu\text{L}$  nuclease-free water to each sample well.
- 13 Seal the wells, then vortex the plate or strip tube for 5 seconds. Verify that all beads have been resuspended, with no visible clumps in the suspension or bead pellets retained on the sides of the wells. Briefly spin to collect the liquid.
- 14 Incubate for 2 minutes at room temperature.
- 15 Put the plate or strip tube in the magnetic stand and leave for 2 minutes or until the solution is clear.

## Step 13. Remove adaptor-dimers with AMPure XP beads (optional)

**16** Remove 25  $\mu\text{L}$  of cleared supernatant to a fresh well. You can discard the beads at this time.

**Stopping Point** If you do not continue to the next step, seal the wells and store at  $-20^{\circ}\text{C}$ .

## 2 Sample Preparation

### Step 14. Prepare samples for multiplexed sequencing

#### Step 14. Prepare samples for multiplexed sequencing

The number of indexed libraries that may be multiplexed in a single sequencing lane is determined by the output specifications of the platform used, together with the amount of sequencing data required for your research design. Calculate the number of indexes that can be combined per lane, according to the capacity of your platform and the amount of sequencing data required per sample.

- 1 Combine the libraries such that each index-tagged sample is present in equimolar amounts in the pool. For each library, use the formula below to determine the amount of indexed sample to use.

$$\text{Volume of Index} = \frac{V(f) \times C(f)}{\# \times C(i)}$$

where  $V(f)$  is the final desired volume of the pool,

$C(f)$  is the desired final concentration of all the DNA in the pool

$\#$  is the number of indexes, and

$C(i)$  is the initial concentration of each indexed sample.

Table 20 shows an example of the amount of 4 index-tagged samples (of different concentrations) and Low TE needed for a final volume of 20  $\mu\text{L}$  at 10 nM.

**Table 20** Example of indexed sample volume calculation for total volume of 20  $\mu\text{L}$

Component	V(f)	C(i)	C(f)	#	Volume to use ( $\mu\text{L}$ )
Sample 1	20 $\mu\text{L}$	20 nM	10 nM	4	2.5
Sample 2	20 $\mu\text{L}$	10 nM	10 nM	4	5
Sample 3	20 $\mu\text{L}$	17 nM	10 nM	4	2.9
Sample 4	20 $\mu\text{L}$	25 nM	10 nM	4	2
Low TE					7.6

- 2 Adjust the final volume of the pooled library to the desired final concentration.

## Step 14. Prepare samples for multiplexed sequencing

- If the final volume of the combined index-tagged samples is less than the desired final volume,  $V(f)$ , add Low TE to bring the volume to the desired level.
  - If the final volume of the combined index-tagged samples is greater than the final desired volume,  $V(f)$ , lyophilize and reconstitute to the desired volume.
- 3** If you store the library before sequencing, add Tween 20 to 0.1% v/v and store at  $-20^{\circ}\text{C}$  short term.

## 2 Sample Preparation

### Step 15. Prepare and analyze sequencing samples

## Step 15. Prepare and analyze sequencing samples

Proceed to cluster amplification using the Illumina Paired-End Cluster Generation Kit; refer to the manufacturer's instructions for this step. The optimal seeding concentration for cluster amplification from SureSelect mRNA libraries is approximately 10-12 pM.

#### NOTE

The optimal seeding concentration may vary, depending on the method used for library quantification and fragment size distribution.

This protocol has been validated with 2 x 100-base paired-end reads. However, read length can be adjusted to achieve the desired research goals.

### Sequencing run setup guidelines for 8-bp indexes

Sequencing runs must be set up to perform an 8-nt index read. For the HiSeq platform, use the *Cycles* settings shown in [Table 21](#). Cycle number settings can be specified on the *Run Configuration* screen of the instrument control software interface after choosing *Custom* from the index type selection buttons.

For complete index sequence information, see [Table 26](#) on page 49.

**Table 21** HiSeq platform Run Configuration screen Cycle Number settings\*

Run Segment	Cycle Number
Read 1	100
Index 1 (i7)	9
Index 2 (i5)	0
Read 2	100

\* Settings apply to v3.0 SBS chemistry.


### Sequence analysis guidelines

The SureSelect<sup>XT</sup> RNA sequencing library preparation method preserves RNA strandedness as described here. The first strand of cDNA is the reverse complement of the poly(A) RNA transcript strand. Since the second strand of cDNA is eliminated before PCR, the sequence of read 1, which starts at the P5 end, matches only the first strand of cDNA. Read 2, which starts at the P7 end, matches the second strand of cDNA (the poly(A) RNA transcript strand). When running analysis of this data to determine strandedness, it is important to include this information. For example, when using the Picard tools (<http://picard.sourceforge.net/>) to calculate RNA sequencing metrics, it is important to include the parameter *STRAND\_SPECIFICITY=SECOND\_READ\_TRANSCRIPTION\_STRAND* to correctly calculate the strand specificity metrics.

## **2 Sample Preparation**

### **Step 15. Prepare and analyze sequencing samples**



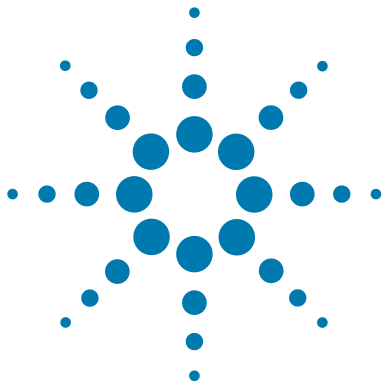

### 3 Reference

Kit Contents 46

Nucleotide Sequences of SureSelect<sup>XT</sup> Indexes A01 to H12 49

This chapter contains reference information, including component kit contents and index sequences.



## Kit Contents

The SureSelect<sup>XT</sup> RNA Reagent Kits contain the following component kits:

**Table 22** SureSelect RNA-Seq Kit Content

Component Kits	Storage Condition	16 Samples	96 Samples
SureSelect Strand Specific RNA Library Prep, ILM, Box 1	-20°C	5500-0134	5500-0135
SureSelect Strand Specific RNA Library Prep, ILM, Box 2	4°C	5190-6410	5190-6411

The contents of each of the component kits listed in [Table 22](#) are described in the tables below.

**Table 23** SureSelect Strand Specific RNA Library Prep, ILM, Box 1 Content

Kit Component	16 Reactions (p/n 5500-0134)	96 Reactions (p/n 5500-0135)
RNA Seq Fragmentation Mix	tube with red cap	bottle
RNA Seq First Strand Master Mix	tube with orange cap	tube with orange cap
RNA Seq Second-Strand + End-Repair Enzyme Mix	tube with blue cap	bottle
RNA Seq Second-Strand + End-Repair Oligo Mix	tube with yellow cap	tube with yellow cap
RNA Seq dA Tailing Master Mix	tube with green cap	bottle
SureSelect Ligation Master Mix	tube with purple cap	tube with purple cap
SureSelect Oligo Adaptor Mix	tube with blue cap	tube with blue cap
RNA Seq PCR Master Mix	tube with red cap	bottle
Uracil DNA Glycosylase (UDG)	tube with yellow cap	tube with yellow cap
SureSelect Primer	tube with brown cap	tube with brown cap
RNA Seq ILM Reverse PCR Primer	tube with black cap	tube with black cap
RNA Seq ILM Post-capture PCR Primer*	tube with green cap	tube with green cap
SureSelect <sup>XT</sup> Indexes, 8 bp <sup>†</sup>	SSEL 8bp Indexes A01 through H02, provided in 16 tubes with white caps	SureSelect 8 bp Indexes A01 through H12, provided in blue 96-well plate <sup>‡</sup>

\* The provided SureSelect ILM Post-capture PCR Primer is not used in the workflow described in this manual. This primer is used in the SureSelect RNA Sequencing Target Enrichment workflow detailed in publication part number G9691-90000.

† See [Table 26](#) on page 49 for index sequences.

‡ See [Table 25](#) on page 48 for a plate map. Although the provided plate contains 96 indexing primers, only indexes A01–H06 should be used for the mRNA library preparation workflow. Wells contain enough volume for two mRNA library preparation reactions per index, using the protocol on [page 33](#).

### 3 Reference

#### Kit Contents

**Table 24** SureSelect Strand Specific RNA Library Prep, ILM, Box 2 Content

Kit Component	16 Reactions	96 Reactions
Oligo(dT) Microparticles	tube with brown cap	bottle
RNA Seq Bead Binding Buffer	tube with purple cap	bottle
RNA Seq Bead Washing Buffer	bottle	bottle
RNA Seq Bead Elution Buffer	tube with green cap	bottle
Nuclease Free Water	bottle	bottle

**Table 25** Plate map for SSEL 8bp Indexes A01 through H12 provided in blue plate in Library Prep kit p/n 5500-0135. Use only indexes A01–H06 (Columns 1 to 6) for the mRNA library preparation workflow.

	1	2	3	4	5	6	7	8	9	10	11	12
A	A01	A02	A03	A04	A05	A06	A07	A08	A09	A10	A11	A12
B	B01	B02	B03	B04	B05	B06	B07	B08	B09	B10	B11	B12
C	C01	C02	C03	C04	C05	C06	C07	C08	C09	C10	C11	C12
D	D01	D02	D03	D04	D05	D06	D07	D08	D09	D10	D11	D12
E	E01	E02	E03	E04	E05	E06	E07	E08	E09	E10	E11	E12
F	F01	F02	F03	F04	F05	F06	F07	F08	F09	F10	F11	F12
G	G01	G02	G03	G04	G05	G06	G07	G08	G09	G10	G11	G12
H	H01	H02	H03	H04	H05	H06	H07	H08	H09	H10	H11	H12

## Nucleotide Sequences of SureSelect<sup>XT</sup> Indexes A01 to H12

Each index is 8 nt in length. See [page 42](#) for sequencing run setup requirements for sequencing libraries using 8-bp indexes.

**Table 26 SureSelect RNA Seq Indexes, for indexing primers in white-capped tubes or blue 96-well plates**

Index	Sequence	Index	Sequence	Index	Sequence	Index	Sequence
A01	ATGCCTAA	A04	AACTCACC	A07	ACGTATCA	A10	AATGTTGC
B01	GAATCTGA	B04	GCTAACGA	B07	GTCTGTCA	B10	TGAAGAGA
C01	AACGTGAT	C04	CAGATCTG	C07	CTAAGGTC	C10	AGATCGCA
D01	CACTTCGA	D04	ATCCTGTA	D07	CGACACAC	D10	AAGAGATC
E01	GCCAAGAC	E04	CTGTAGCC	E07	CCGTGAGA	E10	CAACCACA
F01	GACTAGTA	F04	GCTCGGTA	F07	GTGTTCTA	F10	TGGAACAA
G01	ATTGGCTC	G04	ACACGACC	G07	CAATGGAA	G10	CCTCTATC
H01	GATGAATC	H04	AGTCACTA	H07	AGCACCTC	H10	ACAGATTC
A02	AGCAGGAA	A05	AACGCTTA	A08	CAGCGTTA	A11	CCAGTTCA
B02	GAGCTGAA	B05	GGAGAACA	B08	TAGGATGA	B11	TGGCTTCA
C02	AAACATCG	C05	CATCAAGT	C08	AGTGGTCA	C11	CGACTGGA
D02	GAGTTAGC	D05	AAGGTACA	D08	ACAGCAGA	D11	CAAGACTA
E02	CGAACTTA	E05	CGCTGATC	E08	CATACCAA	E11	CCTCCTGA
F02	GATAGACA	F05	GGTGCGAA	F08	TATCAGCA	F11	TGGTGGTA
G02	AAGGACAC	G05	CCTAATCC	G08	ATAGCGAC	G11	AACAACCA
H02	GACAGTGC	H05	CTGAGCCA	H08	ACGCTCGA	H11	AATCCGTC
A03	ATCATTCC	A06	AGCCATGC	A09	CTCAATGA	A12	CAAGGAGC
B03	GCCACATA	B06	GTACGCAA	B09	TCCGTCTA	B12	TTCACGCA
C03	ACCACTGT	C06	AGTACAAG	C09	AGGCTAAC	C12	CACCTTAC
D03	CTGGCATA	D06	ACATTGGC	D09	CCATCCTC	D12	AAGACGGA
E03	ACCTCCAA	E06	ATTGAGGA	E09	AGATGTAC	E12	ACACAGAA
F03	GCGAGTAA	F06	GTCGTAGA	F09	TCTTCACA	F12	GAACAGGC
G03	ACTATGCA	G06	AGAGTCAA	G09	CCGAAGTA	G12	AACCGAGA
H03	CGGATTGC	H06	CCGACAAC	H09	CGCATACA	H12	ACAAGCTA

[www.agilent.com](http://www.agilent.com)

## In This Book

This guide contains information to run the SureSelect Strand-Specific RNA Library Prep for Illumina Multiplexed Sequencing protocol.

© Agilent Technologies, Inc. 2013-2015,  
2017

Version E0, March 2017



p/n G9691-90010



**Agilent Technologies**